1. 双击打开 ToupView 摄像头软件;



2. 按照图中所示,选择设置单位 (um),选择目前正在使用的显微镜物镜倍数;

🖲 Τοι	upView	200							
File	Edit	View	Browse	Setup	Capture	Image	Proce	ss Layer	Measurem
1			Microme	ter (µm)	▼ 10X	•	100%	•	🖽 🤭 \varkappa
Came	ra		Pixel Nanomet	er (nm)	-	џ 🖂			
	amera L	ist	Microme	ter (µm)		\$			
-			Centimet	r (mm) er (cm)		_			
UC	:MOS03:	100KPA	Meter (m Inch (in))		_			
in Cz	opture 8	v Reso	lution			*			
	ō.	Snap			Record				
	_					_			
Live:						<u> </u>			
Snap):					<u> </u>			
Form	nat:					-			
E Fx	oosure	& Gai	n			¥			
💿 To	oupVie	w							
Fil	e Edit	t Vie	w Brow	vse Se	tup Cap	oture I	mage	Process	Layer M
6		2 🗁	Micro	meter (J	um) 💌	10X	- 1	00%	💽 🖽 🤄
Cam	era					IX			
	Camora	Liet				20X			
	camera	LIST				10X VA	- 11		
U	CMOS)3100K	PA						
_		_	_	_			- 1		
<u>©</u>	Capture	e & Re	solution			^			
		Car			Dee				
4	0	SNa	ab.		Rec	oid			
1									
Live	e:					-			

3.选择直线测量工具;

🐌 ToupView				
File Edit View Browse Setup	Capture Image Process	Layer Measurements Op	itions Window Help	
📴 🛃 🎽 🗁 🛛 Micrometer (µm)	▼ 10X ▼ 40%	💽 🖽 🖑 🗼 🖌 🔸	• 🛛 // 1// 1 • 🗆 🔿 🔿	• @ & • \ I
Camera	🔻 🕂 🖬 🛛 🗸 🗸 🗸	deo [UCMOS03100KPA]		
👩 Camera List	*	0 10	0 200	300
UCMOS03100KPA				
Capture & Resolution	*			
Snap	Record			

在摄像头实时 LIVE 模式下,单击起点位置开始测量,单击终点结束,显示的数值即为实际 距离;(图中测微尺一大格为 100um)



图为测量传感器尖端 LIX 长度;

upView	Browne Setup Capture	Image Process Laver	Measurements Ontions Window	Help		
	Micrometer (um) 🔽 10X	▼ 40% ▼ E		1.000.000000000000000000000000000000000	1 🍐 🗚 🖸 😘 Ŧ 👘	
urement	•	# 10 4 Video [UC]	MO503100KPA]			4
ppearance		0.	100	200 300	400 500	600
me bel Type	 Length					
Width	0					
Color Style	f0000	_				
w1		-				
ulation						
	(333.33, 216.43)	•				
gth	51.52					
e e	-0.03 178.24					
ordinates						
.X .Y	1055					
X	1218	6				
.1	630					
		-				
		-				
		20				
		•		G	- 51. 32μm	
		we have a second				
						1.000
		-				
		400				1000
						1.000
		-				1.
		5				
		8				
_						
asurement	🧕 Camera					
全测	晶线段方 》	上. 选择档:	式下选择此线目	马 按 Delete 键	选择 Ves 即可删图	诠
11/21/12/1	主动权///1	公, 201千庆.	IN I REJERILIENT			21.0
Edit View	Browce Setup Canture	Image Process Laver	Measurements Ontions Window	Help		
	Micrometer (um) 💌 10X	▼ 40% ▼ 1		1.000.0000.00	1 🔬 📶 🚳 🗮 🐨	
ment	•	4 Video [UC	10503100KPA			
earance		0	100	200 300	400 500	600
e June						
Width	0					
Color	ff0000					
1						
2 ulation	+					
	(333.33, 216.43)	•				
th	(384.83, 218.01) 51.52	-				
	-0.03					
dinates	178.24	-				
x	1055	-				
r (685	5				
·	690			~		
			ToupView			
		-				
			Are you sure to r	emove all the selected items?		
		208			- 51- 52um	
				<u>res</u> <u>N</u> o		
		-				
		8				1.000
		-				
		-				
		8				
		-				
		50				
		ō.				